E-Prime
Extensions for
tobiipro

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Psychology Software Tools, Inc.
311 23rd Street Extension, Suite 200
Sharpsburg, PA 15215-2821
Phone 412.449.0078
Fax: 412-449-0079
Email: info@pstnet.com

E-Prime Extensions for Tobii Pro 3.2用户手册
PST-101744
Rev 6

Copyright

Copyright 2019 Psychology Software Tools, Inc. All Rights Reserved.

The information in this document is subject to change without notice. Except as permitted under the United States Copyright Act of 1976, no part of this publication may be reproduced, or distributed in any form or by any means, or stored in a database or retrieval system, without prior written permission of Psychology Software Tools, Inc.

Psychology Software Tools, Inc.
311 23rd Street Extension, Suite 200
Sharpsburg, PA 15215-2821
Phone: 412-449-0078
Fax: 412-449-0079
E-mail: info@pstnet.com
Web: www.pstnet.com

For questions or comments regarding this manual or installation assistance:
Please e-mail us at support@pstnet.com or visit us at <https://support.pstnet.com>

Software Notice: The enclosed software is provided for use by a single user who has purchased the software. The software MAY NOT be reproduced or distributed to others. Unauthorized reproduction and/or sales of the enclosed software may result in criminal and civil prosecution (17 USC 506).

Trademark

Psychology Software Tools, Inc., the Psychology Software Tools, Inc. logo, E-Prime®, E-Prime® logo, images are trademarks or registered trademarks of Psychology Software Tools, Inc. Tobii, Tobii Pro, Tobii Pro Lab, Tobii Pro Studio, and the Tobii Pro logo are trademarks or registered trademarks of Tobii Technology AB. Windows® and Excel® are registered trademarks of Microsoft Corporation in the United States and other countries.

This manual describes the installation procedure for the E-Prime Extensions for Tobii Pro 3.2. Please review the manual completely and thoroughly before beginning the system installation.

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